Caution: For Laboratory Use. A product for research purposes only.

[¹²⁵I]-SAR¹-ILE⁸-ANGIOTENSIN II

Product Number: NEX248

Sar-Arg-Val-[¹²⁵I]Tyr-Ile-His-Pro-Ile

LOT SPECIFIC INFORMATION

TECHNICAL DATA SHEET

		Package Size Information
CALCULATED AS OF:	8-Sep-2014	Package Size
		as of
LOT NUMBER:	EJ01740	17-Oct-2014
		370 kBq
SPECIFIC ACTIVITY:	81.4 TBq/mmol	10 µCi
	2200 Ci/mmol	1.85 MBq
	68.3 MBq/µg	50 µCi
	1846 µCi/µg	3.70 MBq
		100 uCi

RADIOCHEMICAL PURITY: $\geq 95\%$

MOLECULAR WEIGHT: ~1192

PACKAGING: [¹²⁵I]-Sar¹-Ile⁸-Angiotensin II is lyophilized from 0.04M sodium phosphate buffer, pH 4.6, containing 1M glycine, 0.18M sodium chloride and 0.25% BSA. It is shipped ambient.

STABILITY AND STORAGE: The lyophilized [¹²⁵I]-Sar¹-Ile⁸-Angiotensin II should be stored at 4°C or lower. Following reconstitution with distilled water to a concentration of approximately 50 μ Ci/ml on calibration date, aliquot and store at

-20°C or lower. Under these conditions the product is stable and usable for at least eight weeks after fresh lot date.

SPECIFIC ACTIVITY: The initial specific activity of $[^{125}I]$ -Sar¹-Ile⁸-Angiotensin II is 2200 Ci/mmol (81 TBq/mmol),1846 μ Ci/ μ g (68.3 MBq/ μ g). Preparative HPLC is used to separate unlabeled Sar¹-Ile⁸-Angiotensin II from $[^{125}I]$ -Sar¹-Ile⁸-Angiotensin II. Upon decay, $[^{125}I]$ -Sar¹-Ile⁸-Angiotensin II undergoes decay catastrophe and the specific activity remains constant with time. However, it is not known what molecular fragments are generated from the decay event or what functional activity these fragments may have in different assays. References on ^{125}I decay and decay catastrophe of ^{125}I labeled compounds are available.¹⁻⁵

RADIOCHEMICAL PURITY: Initially greater than 95% radiochemically pure as determined by HPLC.

PREPARATIVE PROCEDURE: Sar¹-Ile⁸-Angiotensin II is radioiodinated with no carrier added ¹²⁵I using a modification of the Hunter and Greenwood method⁶ and purified by reversed phase HPLC.

AVAILABILITY: [¹²⁵I]-Sar¹-Ile⁸-Angiotensin II is routinely available from stock. It is prepared fresh and packaged for shipment on the second Monday of each month. Please inquire for larger package sizes.

APPLICATIONS: [¹²⁵I]-Sar¹-Ile⁸-Angiotensin II is an angiotensin receptor antagonist⁷ which binds to a single class of receptors.^{8,9} [¹²⁵I]-Sar¹-Ile⁸-Angiotensin II is useful for studying AII receptors when an antagonist ligand is preferred, for receptor localization studies using autoradiographic techniques^{10,11} and for drug screening programs.

HAZARD WARNING: This product contains a chemical (s) known to the state of California to cause cancer. This product also contains a component which is harmful by contact, ingestion and inhalation. It is irritating to the eyes, skin and respiratory tract and is toxic.

RADIATION UNSHIELDED: 280mR/hr/mCi at vial surface.

REFERENCES:

- 1 Doyle, V.M., Buhler, F.R. and Burgisser, E., Eur. J. Pharm 99 353 (1984).
- 2 Schmidt, J., J. Biol. Chem. 259 1160 (1984).
- 3 Loring, R.H., Jones, S.W., Matthews-Bellinger, J. and Salpeter, M.M., J. Biol. Chem. 257 1418 (1982).
- 4 Berridge, M.S., Jiang, V.W. and Welch, M.J., Rad. Res. 82 467 (1980).
- 5 Charlton, D.E., Rad. Res. 107 163 (1986).
- 6 Hunter, W.M. and Greenwood F.C., Nature 194 495 (1962).
- 7 Khosla, M.C (1972).
- 8 Speth, R.C., Singh, R., Smeby, R.R., Ferrario, C.M. and Husain, A. Neuroendocrinology 38 387 (1984).
- 9 Husain, A., Bumpus, F.M., Desilva, P., and Speth, R.C. Proc. Natl. Acad. Sci. U.S.A. <u>84</u> 2489 (1987).
- 10 McKinley, M.J., Allen, A., Clevers, J., Denton, D.A., and Mendelsohn, F.A.O. Brain Res. 375 373 (1986).
- 11 Millan, M., Aguilera, G., Wynn, P.C., Mendelsohn, F.A.O. and Catt, K.J. Methods Enz. 124 590 (1986).

IODINE-125 DECAY CHART HALF LIFE=60 days

DAYS		0	2	4	6	8	10	12	14	16	18
	0	1.000	0.977	0.955	0.933	0.912	0.891	0.871	0.851	0.831	0.812
	20	0.794	0.776	0.758	0.741	0.724	0.707	0.691	0.675	0.660	0.645
	40	0.630	0.616	0.602	0.588	0.574	0.561	0.548	0.536	0.524	0.512
	60	0.500	0.489	0.477	0.467	0.456	0.445	0.435	0.425	0.416	0.406
	80	0.397	0.388	0.379	0.370	0.362	0.354	0.345	0.338	0.330	0.322
1	100	0.315	0.308	0.301	0.294	0.287	0.281	0.274	0.268	0.262	0.256
1	120	0.250	0.244	0.239	0.233	0.228	0.223	0.218	0.213	0.208	0.203

Radiations: Gamma 35.5 keV (7%), X-ray K alpha 27 KeV (112%), K beta 31 keV (24%)

To obtain the correct radioactive concentration or amount for a date before the calibration date: divide by the decay factor corresponding to the number of days before the calibration date. To obtain the correct radioactive concentration or amount for a date after the calibration date: multiply by the decay factor corresponding to the number of days after the calibration date.

PerkinElmer, Inc. Boston, MA 02118 USA Phone: (800) 762-4000 or (+1) 203-925-4602 www.perkinelmer.com



For a complete listing of our global offices, visit

www.perkinelmer.com/lasoffices